


### Environmental conditions



Operating temperature	15-45°C (59-113°F)
Humidity	20-80%, non-condensing
Max pressure	7 bars (72 psi)

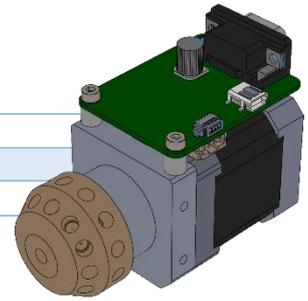

### Fluidic characteristics

Tube port fittings	10-32 UNF, flat bottom
Wetted materials	PTFE (rotor), PEEK (stator)
Channel diameter	0.5 mm (0.020 in) / other size available upon request

### Mechanical characteristics

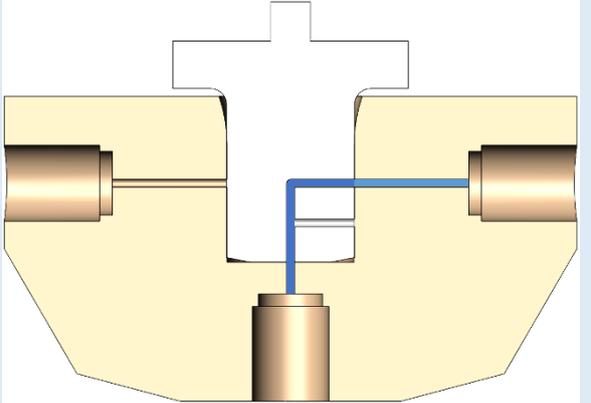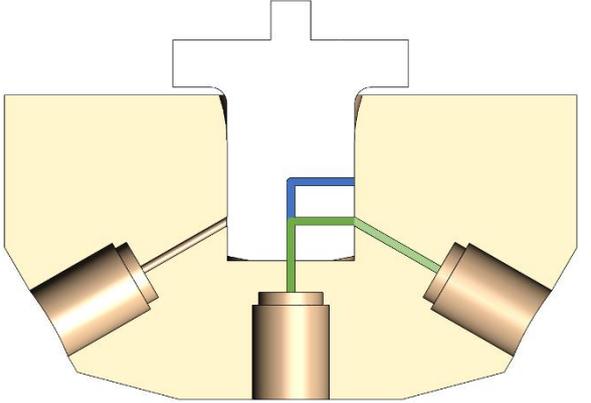Series	LOW CARRYOVER VOLUME
Liquid path (stage 1)	
Liquid path (stage 2)	
<b>Number of channels inside rotary part</b>	2
<b>Number of ports</b>	24
<b>Channel diameter</b>	0.5 mm
<b>Internal volume</b>	3.6 µL
<b>Carryover volume</b>	2.2 µL

**Glossary reminder**


**Internal volume:** Volume inside the system, from entrance to exit.

**Dead volume:** Volume that is “stuck” in the system (dead end), which is not cleanly swept and relies on diffusion to clear out.

**Carryover volume:** Volume of liquid that will be mixed with the next liquid. It is not stuck, but will be swept next time a liquid passes.

COMPLETE MODULE (Ref. P201-O with 24 port valve)

Dimensions	94.8 x 62.8 x 42.3 mm
Weight	0.45 Kg
External fixation system	8 x M3 screws (positions shown in orange below)
Switching time Port to Port	Down to 160 ms

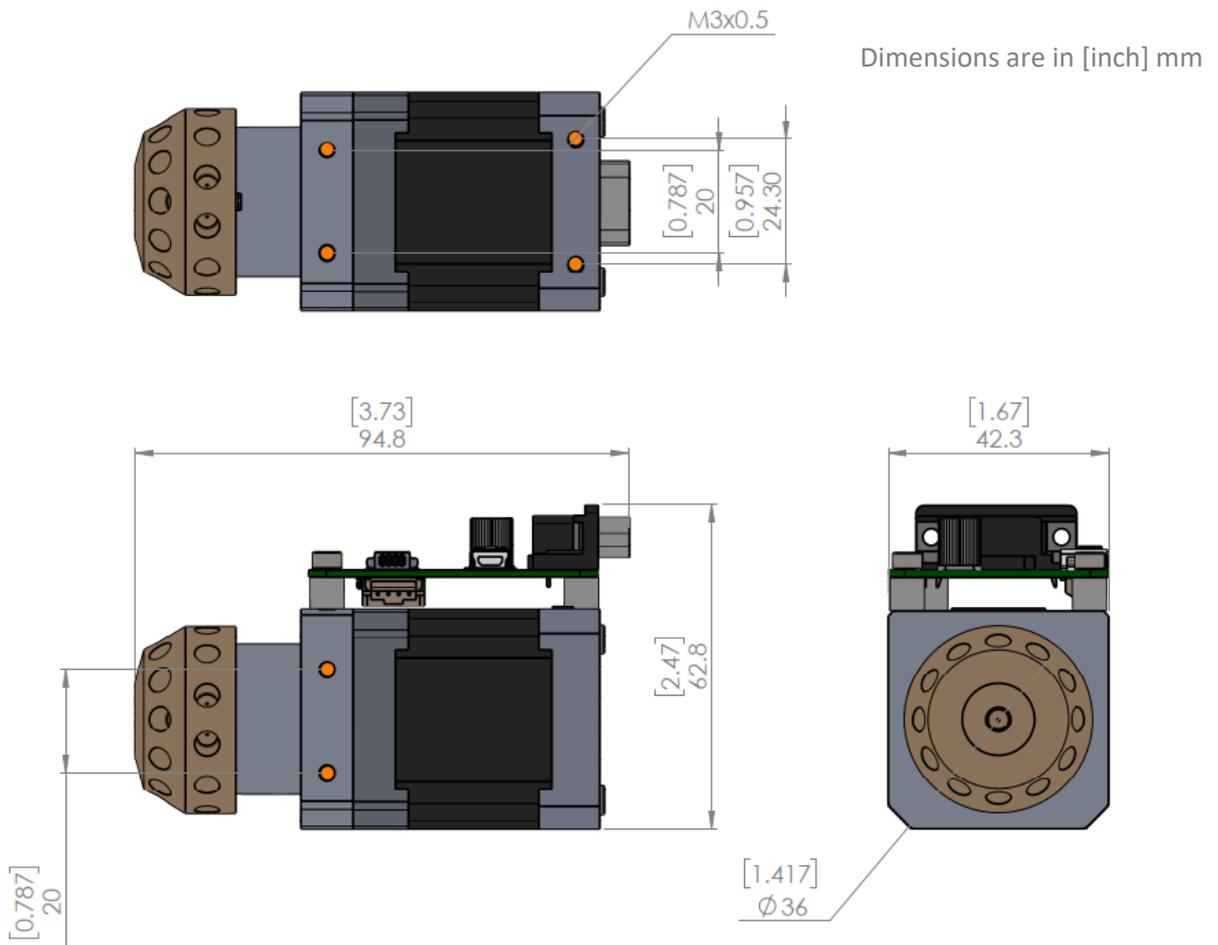

### Electrical characteristics

Power	18-24 VDC, 2.2 A peak
Required cables for operation	USB mini-B male to USB A male or custom cable with PicoBlade™ 4 pos. male, and custom power cord

### Communication and Hardware interface

Interface	USB mini-B, PicoBlade 4 pos., DB9
Communication type	Serial communication on USB mini I2C on PicoBlade™ RS232 on DB9 (RS485 on DB9 upon request) (other upon request)

The valve controller is seen as a virtual serial port. There is an on-board FTDI USB to serial chip, so the drivers are automatically installed on recent Windows 7 and above.

#### SERIAL COMMUNICATION PARAMETERS

Baud rate	9600
Data Bit	8
Parity	None
Stop Bit	1
Handshake	None
End line	<CR>

View the RVM operating manual for the extensive command set.

## Revisions

Version	Author	Release Date	Modifications	Revised By	Approved By
01.01	RRY	04/12/2023	Initial release	MGE	RRY

*The information and data contained in this documentation is subject to change without notice. The reproduction, distribution, and utilization of this document as well as the communication of its contents to others without explicit authorization is prohibited.*

*All rights reserved in the event of the grant of a patent, utility model or design.*

*The general terms and conditions of Advanced Microfluidics SA apply. Alternative agreements must be in written form.*

*Advanced Microfluidics SA shall not be liable for errors contained in this document or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.*